(DE) Gebrauchsinformation

(GB) User instructions

KULZER

MITSUI CHEMICALS GROUP

KULZER GmbH
Leipziger Straße 2
63450 Hanau, Deutschland
Tel. +49 (0) 6181/9689-2574 o. 2571
Fax +49 (0) 6181/9689-3864
technik.wehrheim@kulzer-dental.com
www.kulzer-technik.de

(DE) Gebrauchsinformation Technovit 4002 IQ

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(GB) User Instructions Technovit 4002 IQ

Dear user,

You have received the product Technovit 4002 IQ, a synthetic material that is unique as a two-component, shrinkage-free embedding resin. The product consists of a powder component that is available in your choice of a fast variation (green in colour) or a somewhat slower variation that is white, as well as a slightly viscous liquid.

The materials are subject to the Ordinance on Hazardous Materials. When working with Technovit 4002 IQ, the safety instructions on the packaging must be observed and precautionary measures must be taken that are common when working with chemicals.

Technovit 4002 IQ provides embedding without leaving any gaps and with almost no polymerisation shrinkage. It has outstanding grinding and polishing properties as well as optimum edge definition for the embedded samples.

To ensure these properties and be able to use the embedding system perfectly over its entire service life, it is mandatory to strictly observe several points.

- **Storage of the liquid component at 5 - 21 °C (refrigerator storage recommended)!**

- **Processing of the liquid is recommended in the cooled state!**

- **Avoid direct sunlight – keep in a place that is dark/protected from light!**

Mixing is important for optimum use.

The following parameters are used as a basis for determining the requirements of embedding plastic:

Mixing ratio

1 : 2 to 2 : 3 (Liquid : Powder) by weight

Quantity specification for Kulzer embedding moulds (without sample)

Example embedding mould 40 mm:

1 : 2 = 15 g Liquid : 30 g Powder

2 : 3 = 20 g Liquid : 30 g Powder

If the mixing ratio is too thick, it is possible that bubbles may occur in the mounting resin.

Prepare the liquid and gradually sprinkle in the powder. Thorough mixing is the foundation for optimum embedding results. Mix for 45 seconds, let it stand briefly (swell), mix once again and then pour.

The package sizes are approximately matched to the mixing proportions of the product.

Processing times and curing times (at 21 °C)

Combination of liquid/powder green:
 Processing time 3 min.,
 curing time 10 – 20 min.
 Combination of liquid/powder white:
 Processing time 4 min.,
 curing time 12 – 23 min.

Due to the longer pot life, vacuum infiltration is also possible.

Storage:

Liquid at 5 - 21 °C (refrigerator storage recommended), powder at max. 25 °C.

Delivery units

Technovit 4002 IQ Liquid
 1 x 1,000 ml Art. no. 66064412
 1 x 5,000 ml Art. no. 66064413

Technovit 4002 IQ powder, white
 1 x 1,300 g Art. no. 66064414
 1 x 13,000 g Art. no. 66064415

Technovit 4002 IQ powder, green
 1 x 1,300 g Art. no. 66064416
 1 x 13,000 g Art. no. 66064417

Danger/Safety advice

Please follow the instructions on the product packaging and safety data sheets.

www.kulzer-technik.com

Our technical advice, whether verbal, in writing or by way of trials, is given in good faith but without warranty, and this also applies where proprietary rights of third parties are involved. It does not release you from the obligation to test the products supplied by us as and to their suitability for the intended processes and uses. The application, use and processing of the products are beyond our control, and therefore, entirely your own responsibility. Should, in spite of this, liability be established for any damage, it will be limited to the value of the goods delivered by us and used by you. We will, of course, provide consistent quality of our products within the scope of our General Conditions of Sale and Delivery.

* For more information please visit

www.technovit.com

® = registered trademark,
Kulzer GmbH, Hanau